

## How to install MaxTRAQ 2D with Sentech Camera

Download and install Sentech Camera Drivers

[www.innovision-systems.com/Downloads/Products/Sentech/StCamSWare\\_x64.exe](http://www.innovision-systems.com/Downloads/Products/Sentech/StCamSWare_x64.exe)

Download and install Sentech Camera Support for MaxTRAQ


[www.innovision-systems.com/Downloads/Products/MaxTRAQ/2D/SentechUSB.exe](http://www.innovision-systems.com/Downloads/Products/MaxTRAQ/2D/SentechUSB.exe)

Download and install MaxTRAQ 2D


[www.innovision-systems.com/Downloads/Products/MaxTRAQ/2D/mtsetup.exe](http://www.innovision-systems.com/Downloads/Products/MaxTRAQ/2D/mtsetup.exe)

Once the software has been installed, start MaxTRAQ and select New Recording from the File Menu. This will take you to the recorder view.


1. Select **Sentech USB** as the Camera Type
2. Select the camera you want to use
3. Select Resolution and Frame Rate

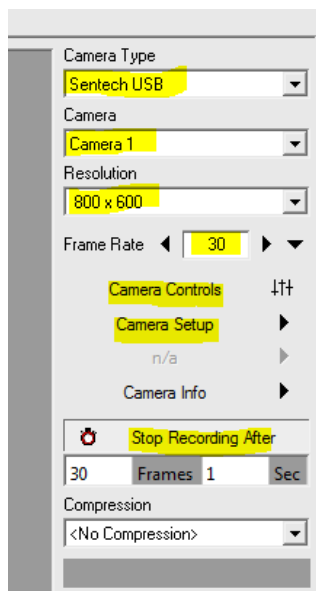

- Click on the viewer button to start live viewing and then click on Camera Controls button and adjust Shutter and Gain

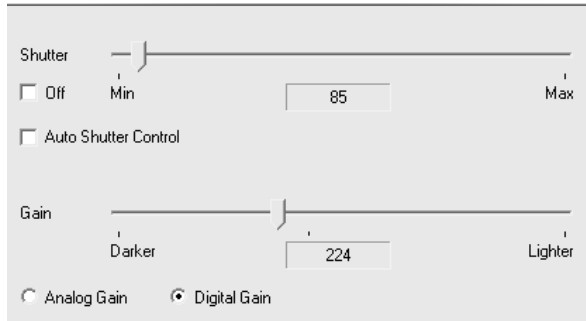

- Click on Camera Setup.

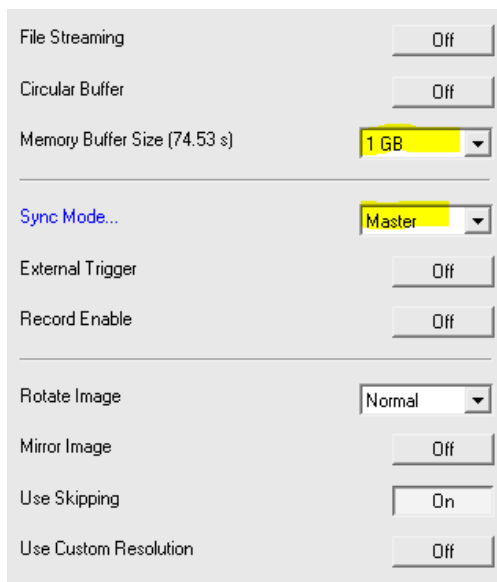

- Select memory buffer size 1 GB
- Sync Mode Master

You can adjust recording time by clicking on Stop Recording After. By default, MaxTRAQ will record until you click stop.

Note; Compression should be set to <No Compression>

To control MaxTRAQ recordings, select Options from the Tools Menu. Please refer to the help for an explanation of the recorder options.