

This document will guide you along in the installation of MaxTRAQ software.

Please Note; If you're planning on using any of our Sentech Cameras with MaxTRAQ, you should install the camera support software prior to installing MaxTRAQ. You can find instructions [here](#).


You must download MaxTRAQ from the Innovision Systems Inc. website at: www.innovision-systems.com.

Once downloaded, double click on the file to launch the installer.


Click **Next** on the first Window.




Choose to accept the license agreement. If you do not agree to the terms of the license agreement, the software will not be installed. Click **Next** after your decision.

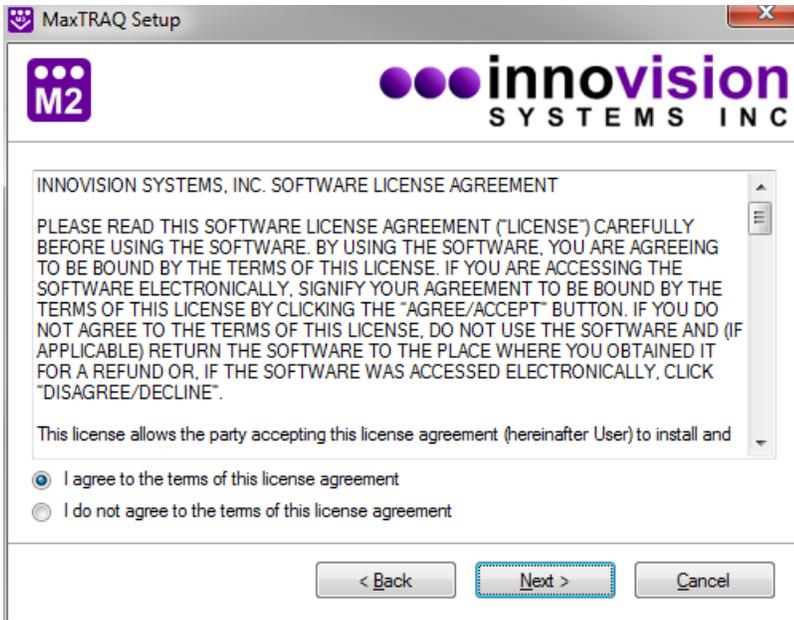


Please do not change the default directories of the installation. This will ease support personnel should you need support. Click **Next**.

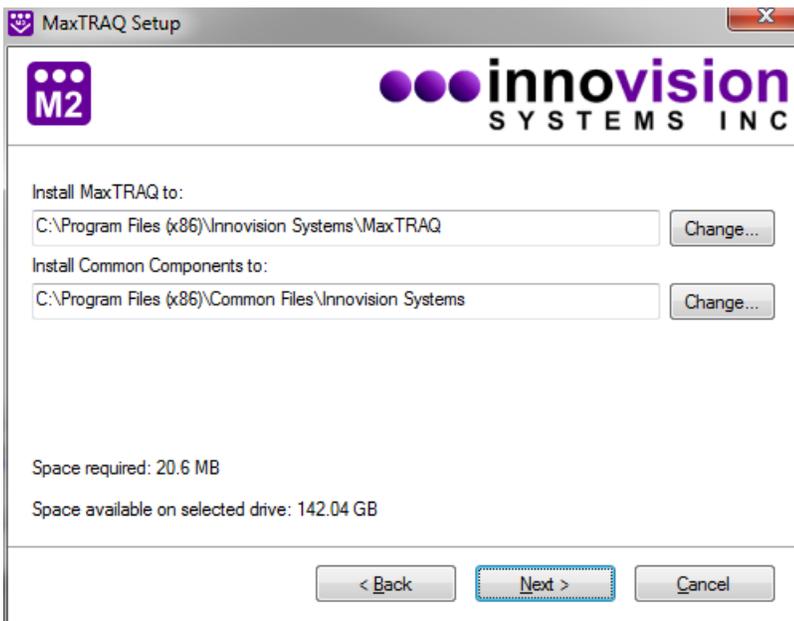


Click on the option you would like for shortcuts and click **Next**.

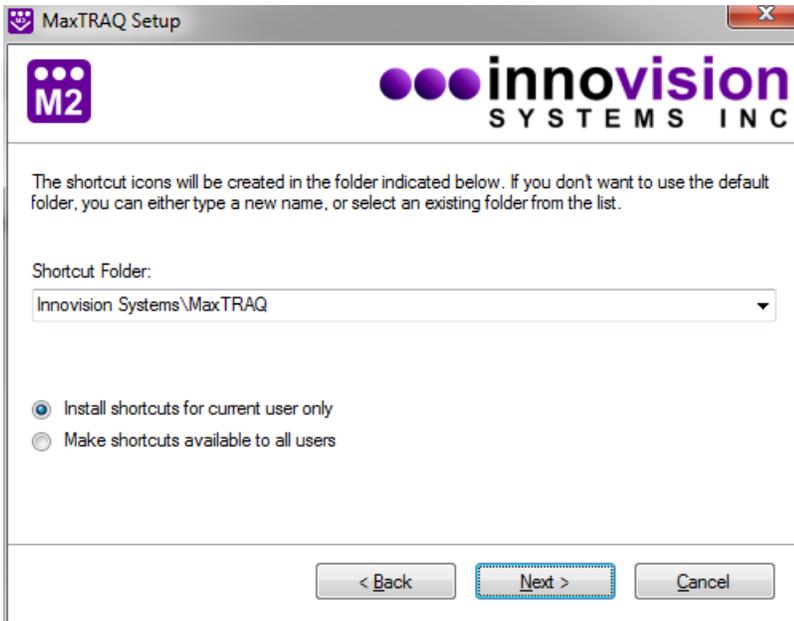

Select one of the installation types. If MaxTRAQ will be used on one computer, select Normal Client Installation. If you are installing on a server to distribute licenses, select Network License Server. Click **Next** after selection. **Note;** You must purchase a network license to ensure the Network License Server functions correctly.

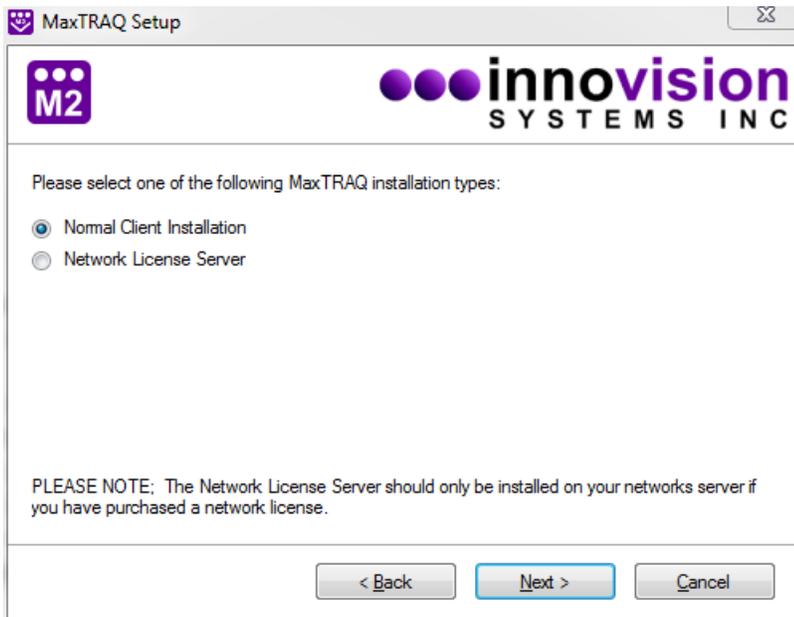


Next you need to select optional components. You can leave the default or if you don't have or plan on using a hardware Dongle for licensing or use MP4 video files you may uncheck either of the options in this dialog. Click the **Next** button.

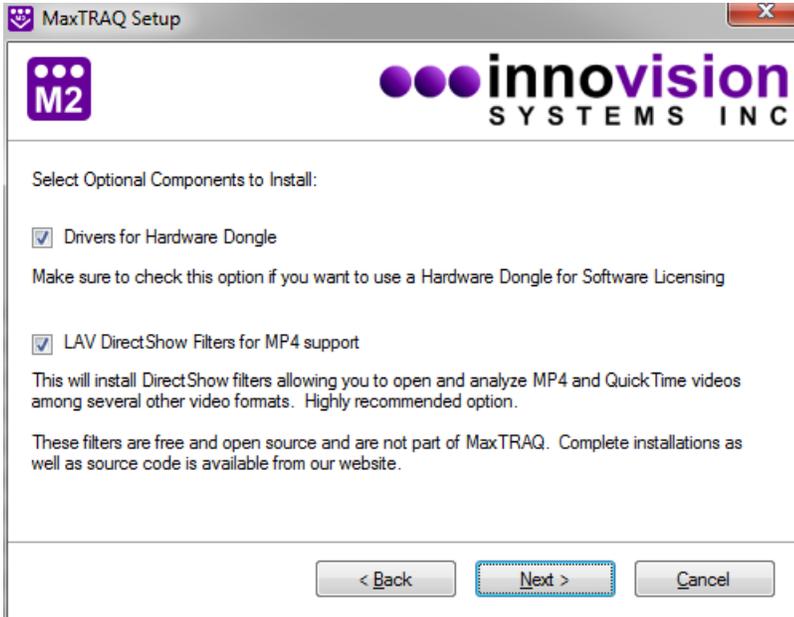

At this window, select the make of camera you will be using, and select **Next**.



Verify that the settings are correct and select **Next**.

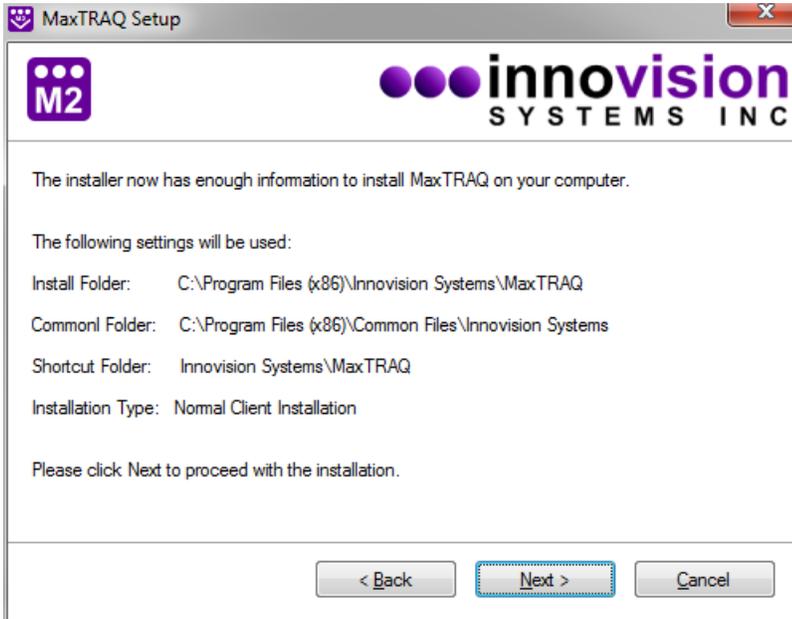

The software install is now complete. Select **Finish** to exit the installer.

